

# NC 40M / 90M

**Laboratory Steam Sterilizers** 





N-s22art



NC 40M / NC 90M laboratory steam sterilizers are specially designed for the steam sterilization of liquid and culture media and also convenient for the sterilization of any labware such as glass and plastics and unpacked solid materials in the laboratories of hospitals, universities and pharmaceutical, food, beverage, dairy and biotechnology industries.

## STANDARD FEATURES FOR LIQUID STERILIZATION

- Flexible temperature sensor to measure the temperature of the liquid load directly as the time required to reach the sterilization temperature of liquid is longer compared to solid loads.
- Cooling function with ambient air ventilation over the chamber that decreases the cooling time of liquid loads to the one third of time spent by conventional systems.
- Condensation of steam at the end of the cycle by means of air ventilated cooling system to prevent a sudden drop in the chamber pressure which causes re-boiling.

#### WELL PRESERVED SENSITIVE MATERIALS

- Second temperature sensor active for solid load sterilization.
- Controlled steam discharge by fractionated exhaust.
- Programmable exhaust rate against sudden temperature and pressure drop to preserve sensitive materials such as glass.







### RELIABLE AND POWERFUL CONTROL SYSTEM: N-Smart™

- Maximum information with highly visible 4,3" colourful LCD display
- Four context-sensitive function keys for versatile & easy use
- Password protected menu to secure run parameters and data
- With 5 Language options: Turkish, English, French, Russian, Spanish
- Icons to make the operations easier for immediate understanding

### **EASY STERILIZATION**

• Five pre-set programs which guarantee safe sterilization of the treated material.


Liquid	121°C	15 minutes
Liquid	121°C	20 minutes
Solid	121°C	20 minutes
Solid	134°C	20 minutes
Solid Quick	134°C	4 minutes

- Two free programs for liquid and solid and a special program for melting purposes which the user can program easily.
- Delayed start function with real time which allows the users to get their samples sterilized on time as they arrive to the laboratory.
- Programmable pre-heating.








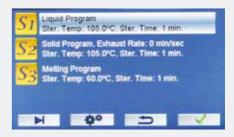


### FULL TRACEABILITY AND REMOTE ACCESS


- Operating parameters can be tracked graphically against time
- Massive data storage with internal memory which stores the records of last 500 cycles detailed and 29.500 cycles as summary
- Memory for the last 100 failures
- USB port to record the operated cycles in the memory to a memory stick
- Possibility of PC Connection through USB / RS 232 port
- Ethernet port for remote access through internet by means of optional NüveCloser™ software
- NuveCloser™ allows access to operating parameters, memory ,failure history and other technical parameters
- Optional panel type printer to keep the records of the sterilization cycle

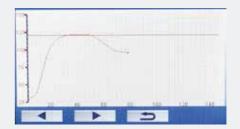






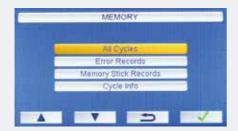

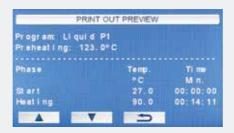

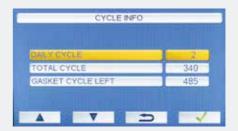



#### SAFETY OVER STANDARDS

- Password protected control system, service and calibration menu
- · Lid safety system
- Lid cannot be opened when the chamber is pressurized
- Program cannot be operated if the lid is open
- Lid can be opened when the liquid temperature is below 80°C after liquid sterilization
- Lid can be opened when the chamber temperature is below 90°C after solid sterilization
- Comprehensive self-diagnostic system to provide the information regarding possible system malfunction. The self-diagnostic system warns user in case of:
- Broken temperature sensors
- Over temperature
- Heater failure
- Low temperature
- Overpressure
- Exhaust tank failure
- Lack of water
- High preheating temperature
- Colour of the display changes to red in case of any alarm condition along with an audible alarm
- Remote failure diagnostic by means of NüveCloser™ software
- Sending e-mails up to five e-mail addresses with the details of failure
- Optional AlerText™ SMS alarm system for ultimate security
- Safety valve against over pressure
- Manometer to follow the chamber pressure
- Chamber manufactured according to PED 2014/68 EC
- Validation port as standard











	NC 40M	NC 90M		
Chamber Volume	40 litres	90 litres		
Temperature Range 105°C - 135°C				
Control System	N-Smart™ Co	ontrol System		
Display	4,3" Colourfu	ıl LCD Display		
No of Preset Programs	2 For Liquids	s, 3 For Solids		
No of Free Programs	1 For Liquids	s, 1 For Solids		
Melting Program Temperature Range	60°C -	100°C		
Timer for Free Programs	1-300 r	ninutes		
Timer for Melting Program	1-60 m	1-60 minutes		
Delayed Start Timer	Programmable With Real Date and Time			
Chamber Material	Stainle	Stainless Steel		
No. of Baskets	2 pcs. (Ø 295x242 mm) 3 pcs. (Ø 360x242 mr			
Power Consumption	2.500 W	7.000 W		
Power Supply	220 V 50/60 Hz.	400 V 50/60 Hz., 3 phases+N+G		
Chamber Dimensions (Diameter x Depth) mm	Ø 320x520	Ø 396x750		
External Dimensions ( WxDxH ) mm.	560x600x985	680x660x1050		
Packing Dimensions ( WxDxH ) mm.	570x660x1200	730x760x1200		
Net / Packed Weight kg.	80 / 91	113 / 134		

#### **Loading Capacity for NC 40M Erlenmeyer Flasks**

252   522		1000	2000	2000	5000 I	
	250 ml	500 ml	1000 ml	2000 ml	3000 ml	5000 ml
	3x8 pcs.	2x4 pcs.	2x3 pcs.	1 pc.	1 pc.	1 pc.

#### **Loading Capacity for NC 40M Schott-Duran Bottles**

250 ml	500 ml	1000 ml	2000 ml	3000 ml	5000 ml
3x12 pcs.	2x8 pcs.	2x5 pcs.	3 pc.	1 pc.	1 pc.

#### **Loading Capacity for NC 90M Erlenmeyer Flasks**

	250 ml	500 ml	1000 ml	2000 ml	3000 ml	5000 ml
Γ	4x12 pcs.	3x7 pcs.	3x4 pcs.	2x3 pcs.	2x1 pc.	1 pc.

#### **Loading Capacity for NC 90M Schott-Duran Bottles**

250 ml	500 ml	1000 ml	2000 ml	3000 ml	5000 ml	10000 ml
4x18 pcs.	3x11 pcs.	3x8 pcs.	2x4 pcs.	2x1 pc.	2x1 pc.	1 pc.

#### **FACTORY FITTED OPTIONS**

**ACCESSORIES FOR OT 40L** NC XXMY Panel mounted thermal printer S 09 028 Basket (Height: 345 mm) NuveLift™ to load/unload basket for NC 90M NC 90ME S 09 027 Basket (Height: 242 mm)

**OPTIONS** 

AlerText™ GSM alarm module A 08 191

A 08 212 NuveCloser<sup>™</sup> Software CD with 3 m. RS 232 cable

#### S 09 040 Basket (Height: 160 mm) (Unperforated)

**ACCESSORIES FOR OT 90L** 

Basket (Height: 360 mm) S 09 034 S 09 035 Basket (Height: 242 mm) S 09 041 Basket (Height: 170 mm)

Basket (Height: 242 mm) (Special for NC 90ME) S 09 043

